



MAINTENANCE Manual

for Laboratory Equipment

2nd Edition





MAINTENANCE Manual

for Laboratory Equipment

2nd Edition

WHO Library Cataloguing-in-Publication Data

Maintenance manual for laboratory equipment, 2nd ed.

1.Laboratory equipment. 2.Maintenance. 3.Manuals. I.World Health Organization. II.Pan American Health Organization.

ISBN 978 92 4 159635 0 (NLM classification: WX 147)

© World Health Organization 2008

All rights reserved. Publications of the World Health Organization can be obtained from WHO Press, World Health Organization, 20 Avenue Appia, 1211 Geneva 27, Switzerland (tel.: +41 22 791 3264; fax: +41 22 791 4857; e-mail: bookorders@who.int). Requests for permission to reproduce or translate WHO publications – whether for sale or for noncommercial distribution – should be addressed to WHO Press, at the above address (fax: +41 22 791 4806; e-mail: permissions@who.int).

The designations employed and the presentation of the material in this publication do not imply the expression of any opinion whatsoever on the part of the World Health Organization concerning the legal status of any country, territory, city or area or of its authorities, or concerning the delimitation of its frontiers or boundaries. Dotted lines on maps represent approximate border lines for which there may not yet be full agreement.

The mention of specific companies or of certain manufacturers' products does not imply that they are endorsed or recommended by the World Health Organization in preference to others of a similar nature that are not mentioned. Errors and omissions excepted, the names of proprietary products are distinguished by initial capital letters.

All reasonable precautions have been taken by the World Health Organization to verify the information contained in this publication. However, the published material is being distributed without warranty of any kind, either expressed or implied. The responsibility for the interpretation and use of the material lies with the reader. In no event shall the World Health Organization be liable for damages arising from its use.

Design and Layout: L'IV Com Sàrl, Morges Switzerland

Printed in Spain

Contact:

Dr G. Vercauteren, Coordinator, Diagnostics and Laboratory Technology, Department of Essential Health Technologies, World Health Organization, 20 Avenue Appia, 1211 Geneva 2, Switzerland

This document is available at www.who.int/diagnostics_laboratory



Table of Contents

TABLE OF FIGURES	viii
ACKNOWLEDGEMENTS	x
INTRODUCTION	xi
CHAPTER 1 • MICROPLATE READER	1
Photograph of microplate reader	1
Purpose of the microplate reader	1
Operation principles	1
Installation requirements	3
Routine maintenance	3
Troubleshooting table	4
Basic definitions	5
CHAPTER 2 • MICROPLATE WASHER	7
Photograph of microplate washer	7
Purpose of the microplate washer	7
Operation principles	7
Installation requirements	9
Routine maintenance	9
Troubleshooting table	11
Basic definitions	12
CHAPTER 3 • pH METER	13
Purpose of the equipment	13
Photograph and components of the pH meter	13
Operation principles	13
pH meter components	14
Typical circuit	15
Installation requirements	16
General calibration procedure	16
General maintenance of the pH meter	17
Basic maintenance of the electrode	18
Troubleshooting table	18
Basic definitions	19
Annex: The pH theory	20

CHAPTER 4 • BALANCES	21
Photographs of balances	21
Purpose of the balance	22
Operation principles	22
Installation requirements	26
Routine maintenance	27
Troubleshooting table	28
Basic definitions	29
 CHAPTER 5 • WATER BATH	 31
Diagram of a water bath	31
Operation principles	31
Water bath controls	32
Water bath operation	32
Troubleshooting table	34
Basic definitions	34
 CHAPTER 6 • BIOLOGICAL SAFETY CABINET	 35
Illustration of a biological safety cabinet	35
Purposes of the equipment	35
Operation principles	35
Biological safety	39
Installation requirements	39
Using the safety cabinet	39
Routine maintenance	40
Functional evaluation (alternative)	41
Table of functional evaluation of biological safety cabinets	42
Troubleshooting table	43
Basic definitions	44
 CHAPTER 7 • CENTRIFUGE	 45
Photograph of centrifuge	45
Purpose of the centrifuge	45
Operation principles	45
Components of the centrifuge	46
Installation requirements	48
Routine maintenance	48
Appropriate management and storage recommendations	48
Troubleshooting table	50
Basic definitions	52
 CHAPTER 8 • WATER DISTILLER	 53
Diagram of a water distiller	53
Purpose of the water distiller	53
Operation principles	54
Installation requirements	54
Routine maintenance	55
Troubleshooting table	56
Basic definitions	57

CHAPTER 9 • DILUTOR	59
Diagram of a dilutor	59
Purpose of the dilutor	59
Operation principles	60
Installation requirements	61
Routine maintenance	61
Troubleshooting table	63
Basic definitions	64
 CHAPTER 10 • DISPENSER	 65
Photograph and diagram of the dispenser	65
Purpose of the dispenser	65
Requirements for operation	67
Routine maintenance	67
Troubleshooting table	68
Basic definitions	68
 CHAPTER 11 • SPECTROPHOTOMETER	 69
Photograph of spectrophotometer	69
Purpose of the equipment	69
Operation principles	69
Spectrophotometer components	72
Installation requirements	73
Spectrophotometer maintenance	73
Good practices when using the spectrophotometer	75
Troubleshooting table	77
Basic definitions	79
 CHAPTER 12 • AUTOCLAVE	 81
Photograph of the autoclave	81
Purpose of the autoclave	81
Operation principles	82
Operation of the autoclave	84
Installation requirements	87
Routine maintenance	88
Maintenance of specialized components	90
Troubleshooting table	91
Basic definitions	92
 CHAPTER 13 • DRYING OVEN	 93
Photograph of drying oven	93
Purpose of the oven	93
Operating principles	93
Installation requirements	94
Oven operation	94
Oven controls	95
Quality control	96
Routine maintenance	96
Troubleshooting table	97
Basic definitions	98

CHAPTER 14 • INCUBATOR	99
Photograph of incubator	99
Operating principles	99
Incubator controls	101
Installation requirements	101
Routine maintenance and use of the incubator	101
Troubleshooting table	103
Basic definitions	104
 CHAPTER 15 • MICROSCOPE	 105
Photographs of microscopes	105
Purpose of the equipment	106
Operation principles	106
Installation requirements	108
Description of potential problems with microscopes	109
General maintenance of the microscope	111
Troubleshooting table	115
Basic definitions	116
 CHAPTER 16 • PIPETTES	 119
Photographs of pipettes	119
Purpose of the pipettes	120
Operation principles of the pipette	120
Requirements for use	120
Using the pipette	121
Routine maintenance	122
Troubleshooting table	125
Basic definitions	126
 CHAPTER 17 • STIRRING HEATING PLATE	 127
Photograph of the stirring heating plate	127
Operation principles	127
Controls of the stirring heating plate	127
Installation requirements	128
Operation of the stirring heating plate	128
Routine maintenance	128
Troubleshooting table	129
Basic definitions	129
 CHAPTER 18 • REFRIGERATORS AND FREEZERS	 131
Photograph of a refrigerated storage unit	131
Purpose of refrigerated storage units	132
Operation principles	132
Installation requirements	133
Refrigerator control circuit	134
Refrigerator operation	134
Refrigerator routine maintenance	135
Troubleshooting table	137

Operation of ultralow freezers	138
Turning the unit on	138
Routine maintenance	139
Troubleshooting table	140
Basic definitions	141

CHAPTER 19 • CHEMISTRY ANALYSERS	143
Photographs of chemistry analysers	143
Purpose of chemistry analysers	144
Operation principle	144
Components	144
Installation requirements	145
Operation of the dry chemistry analyser	145
Operation of the wet chemistry analyser	146
Routine maintenance of chemistry analysers	146
Non-routine maintenance and troubleshooting	147
Troubleshooting table	148
Basic definitions	148

CHAPTER 20 • COLORIMETERS	149
Photograph of colorimeter	149
Purpose of the colorimeter	149
Operating principle	149
Components	150
Installation requirements	150
Operation of the colorimeter	150
Operation of the haemoglobinometer	151
Routine maintenance	151
Troubleshooting table	154
Basic definitions	155

BIBLIOGRAPHY	157
---------------------	-----

预览已结束，完整报告链接和二维码如下：

https://www.yunbaogao.cn/report/index/report?reportId=5_29500

